








IndiMag[®] 2




Safety Sheet




WARNING	
	Bypassing or putting safety devices out of operation is strictly prohibited and constitutes improper use.
	Housing parts and covers must be closed during operation and may only be opened to eliminate malfunctions.
	Danger of crushing! Reaching into the unit during operation can lead to injuries.


WARNING	
	Electric shock hazard! Maintenance and repair work may only be carried out when the unit is disconnected from the power supply and the front power switch stops glowing.


WARNING	
	Warning: Biological hazard Use caution when loading or removing infectious samples into or from the IndiMag 2. Handle biological samples as if they may contain potentially infectious agents.
	Warning: Risk of contamination Contain and clean contamination from spilled samples immediately.


WARNING	
	Warning! Hazardous chemicals! Do not use bleach to wipe the device on the inside.


NOTICE	
	Avoid spraying cleaning agents directly into the device.
	Do not use acetone to wipe the front screen.
	Caution: Risk of damage to the instrument Avoid spilling chemicals or other liquids onto or into the IndiMag 2. Damage caused by liquid spillage will void the warranty.

WARNING	
	Danger of crushing when transporting the unit.
	If the unit is not transported properly, this can lead to personal and serious injury. <ul style="list-style-type: none"> • Only move the unit in an upright position. Do not tilt the unit. • Only lift the unit as far off the ground as necessary. • Wear personal protective equipment.
	Use personal protective equipment (safety shoes, gloves).

WARNING	
	Burn hazard: Hot surface! Even after a process run the lysis and elution blocks can be hot. Handle with care.

WARNING	
	Warning: Optical radiation! Do not damage the front screen or remove any body sheets. Do not under any circumstances disable the magnetic hood sensors.

NOTICE	
	Ensure that all installation and supply connections of the unit are freely accessible after installation. Ensure that space is left on both sides and behind the IndiMag 2 to allow adequate air circulation.

NOTICE	
	Only original spare parts must be used.



Further information and support

- IndiMag 2 Handbook: www.indical.com/indimag2
- Technical assistance: support@indical.com

For up-to-date licensing information and product-specific disclaimers, see the respective INDICAL kit handbook or user manual. Trademarks: IndiMag[®] and INDICAL[®] (INDICAL BIOSCIENCE). Registered names, trademarks, etc., used in this document, even when not specifically marked as such, are not to be considered unprotected by law. HB-2662-EN-002 05/2024 © 2024 INDICAL BIOSCIENCE GmbH, all rights reserved.