

IndiMag® Handheld Quick Start Protocol


For manual processing of magnetic bead-based extraction kits, e.g. the IndiMag Pathogen IM48 Cartridge (SP947654P608) or the IndiMag Pathogen Kit w/o plastics (cat. no. SP947257)

Further information and support


- IndiMag Pathogen Kit w/o plastics, IndiMag Pathogen IM48 Cartridge and IndiMag Handheld Handbook: www.indical.com/handbooks
- Technical assistance: support@indical.com

For up-to-date licensing information and product-specific disclaimers, see the respective INDICAL kit handbook or user manual. Trademarks: IndiMag® and INDICAL® (INDICAL BIOSCIENCE); Torx® (Acument Intellectual Properties, LLC, MI, USA). Registered names, trademarks, etc., used in this document, even when not specifically marked as such, are not to be considered unprotected by law. HB-2702-EN-001 12/2025 © 2025 INDICAL BIOSCIENCE GmbH, all rights reserved.

Assembly

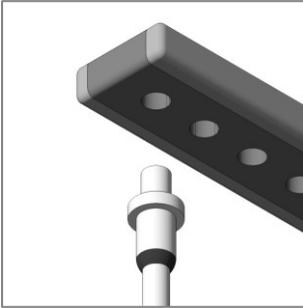

1. Insert the thicker end of the magnet rod into the opening of the IndiMag Handheld.

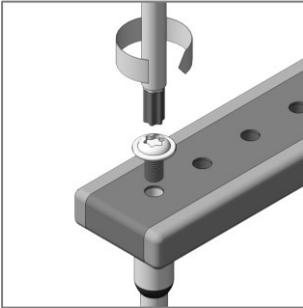

2. Place screw through IndiMag Handheld and tighten it using the T10 Torx® angle key wrench.
3. Repeat the procedure until all eight magnetic rods are incorporated into the IndiMag Handheld.

Important: After assembly, check and ensure all eight magnetic rods are levelled by placing the rod-end on a flat surface (e.g., workbench) and check if all rods are touching the flat surface.

Note: Always place a rod cover (IndiMag 48 PW Rod cover, cat. no. PW940237) when using the IndiMag Handheld so it covers all eight magnetic rods.

CAUTION! The IndiMag Handheld contains strong permanent magnets. Keep loose ferrous objects away and avoid direct contact with the magnet material.

CAUTION! Persons with pacemakers or other implanted medical devices must keep a safe distance, as magnetic fields may interfere with their function.

CAUTION! Keep the device away from magnet-sensitive items such as watches, credit cards, magnetic tapes, and electronic equipment. Maintain distance between multiple units.

Change index

PCard	Version	Change
HB-2702-EN-001	December 2025	Product launch