

Prefilled IndiMag Pathogen KF96 Cartridges

Ready-to-load reagent plates for automated DNA/RNA extraction, compatible with KingFisher™ and equivalent 96-well platforms.

Peel. Add sample & lysis buffer. Run.

Prefilled, sealed, and ready to run

- Preloaded beads and buffers

Increase workflow efficiency

- Skip the pipetting marathon
- Reduced risk of errors with fewer steps

Performance without the plastic burden

- No need to order or store extra plastics
- Reduced reagent bottle waste




Get in touch for more information

Support: us_support@indical.com | Orders: us_orders@indical.com | Call: +1 (407) 766 5128


Simplified DNA/RNA Extraction with IndiMag Pathogen KF96 Cartridges



Shake, spin, and peel off the foil

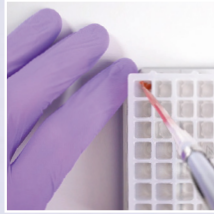

Add sample and Lysis Buffer

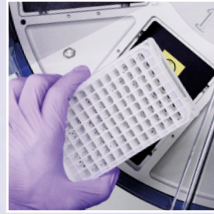

Load and go

IndiMag prefilled cartridges are ready-to-use reagent plates for DNA/RNA extraction, containing magnetic beads, wash and elution buffers sealed in individual wells. No reagent preparation. No additional plastics.

What's in the box?



- A Lysis plate**
Pre-loaded with Proteinase K.
Add sample and Lysis Buffer
- B Wash 1 plate**
Pre-loaded with magnetic beads and wash buffer
- C Wash 2 plate**
Pre-loaded wash buffer
- D Wash 3 plate**
Pre-loaded wash buffer
- E Lysis/binding buffer bottle**
Contains lysis buffer.
Add to lysis plate along with sample.
- F Rod Cover plate**
Ready to load
- G Elution plate**
Pre-loaded with elution buffer; purified DNA/RNA is collected here

Specification	Details
Compatible Instruments	KingFisher Flex, KingFisher Apex, AutoPure 96, BioSprint 96, and equivalent 96-well magnetic particle processors
Sample Types	Whole blood, serum, plasma, swabs, washes, feces, oral/body fluids, tissue and others* <i>*Pretreatment protocols for additional sample types are available by contacting us_support@indical.com.</i>
Target	Viral DNA/RNA and bacterial DNA
Storage & Shipping	Room temperature (15–25 °C / 59–77 °F)

Designed to grow with your lab. Start small, scale fast with flexible kit sizes tailored to your throughput, from 1 x 96 preps to bulk supply.

Product	Cat. no. 96 Preps	Cat. no. 384 Preps	Cat. no. 1536 Preps
IndiMag Pathogen KF96 Cartridge	SP947855P196 (1 x 96)	SP947855P496 (4 x 96)	SP947855P1696 (16 x 96)



Request a quote online: shop.indical.com

Email: us_support@indical.com | us_orders@indical.com | Call: +1 (407) 766 5128

For up-to-date licensing information and product-specific disclaimers, see the respective handbook or user manual. Regulatory requirements vary by country, products may not be available in your geographic area. Trademarks: INDICAL®, IndiMag®, IndiSpin® (INDICAL BIOSCIENCE GmbH). KingFisher™ (Thermo Fisher Scientific). Registered names, trademarks, etc., used in this document, even when not specifically marked as such, are not to be considered unprotected by law. INFLY-IndiMag-KF96-EN-NA-260630 © 2026 INDICAL BIOSCIENCE

INDICAL
BIOSCIENCE

INDICAL Inc.
1317 Edgewater Dr #3722
Orlando, FL 32804, USA